

A Reliable Research Partner in Life Science and Medicine

# **Annexin V-FITC Azid-Free Lyophilized Powder**

Cat. No: E-CK-A111U Size: 25 µg/50 µg/100 µg/200 µg

Cat.	Products	25 μg	50 μg	100 µg	200 μg	Storage
E-CK-A111U	Annexin V-FITC Azide-Free	25 μg	25 μg×2	100 μg	100 μg×2	-20 °C, shading light
	Lyophilized Powder					


## **Storage**

The lyophilized powder is stable for at least two years when stored at  $-20 \, \mathbb{C}$  in the dark after arrival.

#### Introduction

Elabscience® Annexin V-FITC Azide-Free Lyophilized Powder is developed for long-term storage and userdefined concentrations of FITC-conjugated Annexin V. It contains no preservatives like azides or Proclin. Annexin V is a member of the annexin family, which binds to phosphatidylserine (PS) in a calcium-dependent manner. The FITC-conjugated format of this protein, Annexin V-FITC, binds specifically to the PS displayed on the outer leaflet of apoptotic cell membrane. The signal of Annexin V-FITC can be collected with flow cytometry or fluorescence microscopy to detect cell apoptosis.

Jurkat cells were treated with 5 μM Camptothecin and detected with this reagent and PI.

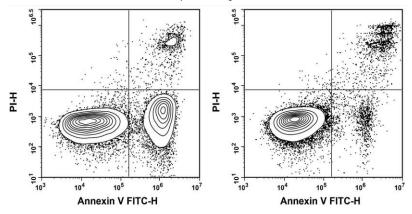

Jurkat cells were cultured with (**Left**) or without (**Right**) 5 μM Camptothecin for 4 h. Annexin V-FITC single-positive cells were early apoptotic cells, Annexin V-FITC and PI double-positive cells were necrotic or late apoptotic cells, and PI single-positive cells were naked nuclei.

#### Reconstitution

Reconstitute the reagent with sterile deionized water. Make sure the final concentration is less than 1 mg/mL. (This concentration is high enough for stock solution. In our internal tests, working solutions with  $1\sim2~\mu g/mL$  Annexin V-FITC have best results for flow cytometry.) For example, at least  $10~\mu L$  water should be added to  $10~\mu g$  Annexin V-FITC. Otherwise, the lyophilized cake may not be thoroughly soaked in water.

### **Cautions**

- 1. This kit is for reseach use only.
- 2. Detect apoptosis as soon as possible after staining to avoid the increase number of apoptosis or necrosis. Avoid extended exposure of the samples to direct light to protect the fluorophores from quenching.
- 3. For your safety and health, please wear the lab coat and disposable gloves before the experiments.

#### For Research Use Only

Toll-free: 1-888-852-8623 Tel: 1-832-243-6086 Fax: 1-832-243-6017

Web: <a href="mailto:www.elabscience.com">www.elabscience.com</a>
<a href="mailto:techsupport@elabscience.com">Email: techsupport@elabscience.com</a>